



Purpose & Scope

The purpose of this document is to provide business advantages, system administrator installation, and end-user access procedures for the use of Remote Desktop (RD) Web Access for Instrument Manager (IM).

Contents

Overview of RD Web Access.....	2
What is RD Web Access?	2
What are the benefits of RD Web Access versus thin client?.....	2
Before You Begin	3
Understanding remote vs. thin client technologies	3
Verifying system requirements	4
Server	4
Client PC.....	4
Setting up RD Web Access for IM	5
Overview of access methods	5
Overview of Windows Installation access	5
Overview of Web URL access.....	5
Set up the Server	6
To install RD Web Access:.....	6
To enable RD Web Access:	6
Set up the Windows Install:	7
To create a Windows Install .msi file:.....	7
To install the Windows Install .msi file on the client PC:.....	7
To verify the Windows Install setup:.....	8
Set up the Web URL	8
To create an .rdp file:	8
To enable the client PC to connect to the server:.....	9
To verify the Web URL setup:	9
Accessing IM with RD Web Access	10
Windows Install	10
To access IM with RD Web Access via a Windows Install:.....	10
Web URL.....	10
To access IM with RD Web Access via a Web URL:	10



Overview of RD Web Access

What is RD Web Access?

Remote Desktop Web Access (RD Web Access) is a technology from Microsoft that provides an alternative deployment method for users to access Instrument Manager (IM) other than the traditional IM thin client method. It uses Microsoft Remote Desktop Web Connection to provide user access to IM via a web browser.

This Microsoft technology allows users to access an application installed on another PC through an internet connection. Remote Desktop Web Connection, available when using RD Web Access, allows users to access an application running on another PC with appropriate permissions.

What are the benefits of RD Web Access versus traditional thin client installation?

Benefits	Details
Faster Deployment Time and Reduced Maintenance	Whether it is for initial deployment or upgrading to a new version, it eliminates the need to install or upgrade IM thin client software on each computer system that requires access to IM.
Streamlined Installation	Unlike IM thin client installations, remote desktop technology requires neither modification of the IMOPTION.ini file, nor sharing of an IM directory on the server.
Increased Deployment Scalability	<p>You can use remote desktop technology to eliminate the need to update numerous IM thin client PCs manually during installation or upgrades:</p> <ul style="list-style-type: none"> • You can connect to IM using only your Internet Explorer browser on a client PC. • Or, you can create a Windows install file, which is typically much smaller (< 700 KB) than a full thin client IM installation. Using remote desktop technology, the software can be pushed out and run on a client PC as part of your organization's IT policy.
Secured Connectivity and Reduced Bandwidth Requirements	<p>Data can be optionally secured via SSL certificate. Further, data security is improved because data is never physically sent to or stored on the client PC. Instead of transferring data to thin client PCs for display, network traffic with remote desktop technology consists only of "screen scrape" data—so data is live on the server PC only.</p> <p>Better security as File and Print Sharing do not have to be enabled on the server to access Instrument Manager.</p>



Before You Begin


Understanding remote desktop vs. thin client technologies

Before you set up remote desktop technology, it is recommended that you review IM thin client concepts, installation, system requirements, and administrator workflows, which are described in the *Instrument Manager Installer's Guide*.

In addition, for an overview about remote desktop technologies, see Microsoft's website:

<http://technet.microsoft.com/en-us/library/dd759202.aspx>

This high-level network diagram compares instances of full IM installations, depending upon thin client or remote technologies:

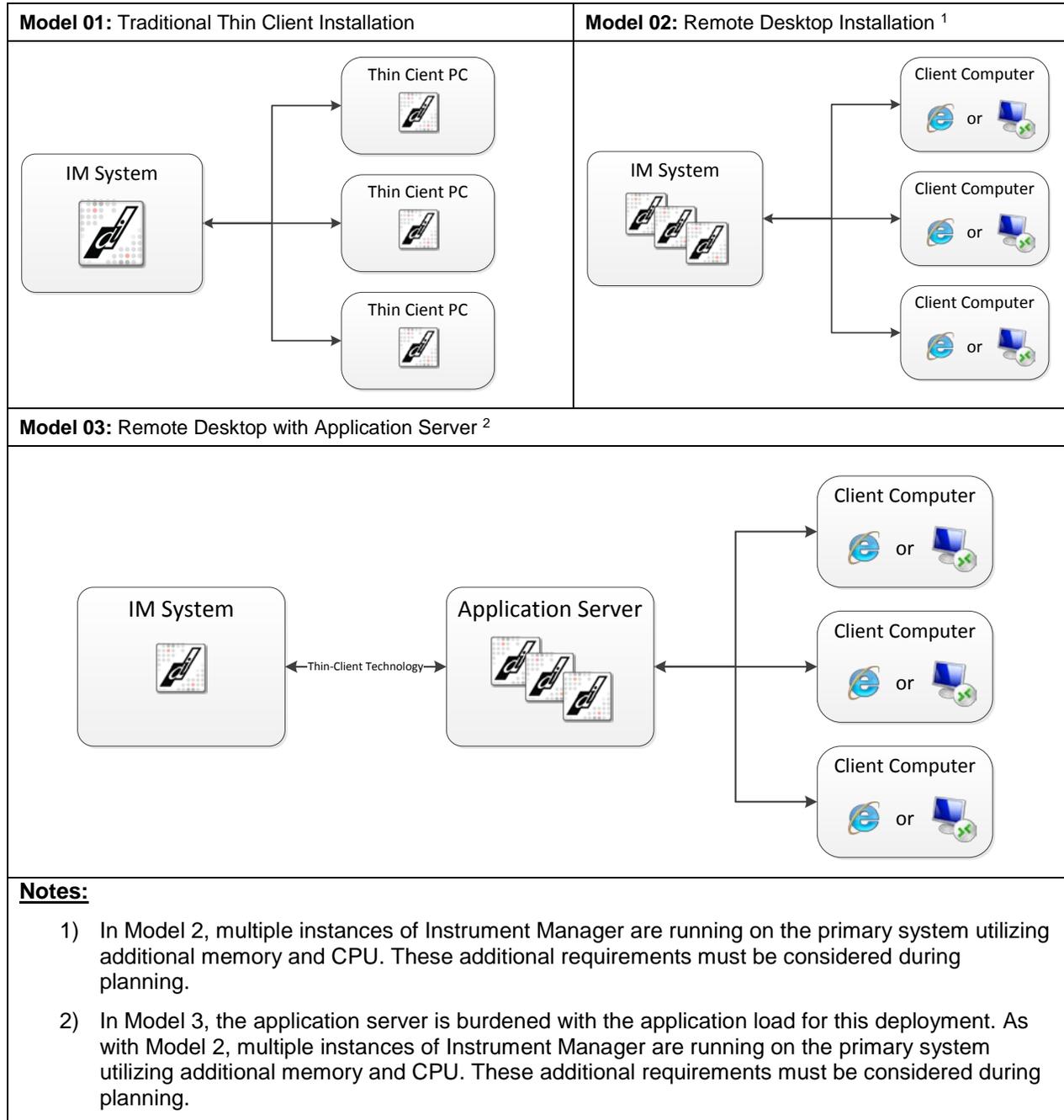



Verifying System Requirements

Primary System or Server

In order to use RD Web Access for IM, the primary IM system must have the following requirements:

- Windows 2008 R2 Server (64-bit) or a Windows 7 (64-bit) system
- A digital security certificate if you wish to apply it to the .rdp file for web URL access
- 256MB of additional Memory needed per client PC connected (depends on functionality used within Instrument Manager)

The following is needed to perform RD Web Access setup:

- Server Login Credentials (Windows login and password)
- Server Host Name

Client PC

Confirm system requirements for the client PC that will use RD Web Access:

- Microsoft Internet Explorer v6.0 or greater is required for best ActiveX Client Control performance.
Note: Other browsers such as Google Chrome, Mozilla Firefox, etc... maybe compatible with Microsoft's Remote Desktop Web (RD Web) but have not been tested and capabilities and/or restrictions are unknown.
- Client computers must be have a version f Remote Desktop Connection (RDC) that supports at least Remote Desktop Protocol (RDP) 6.1.
- ActiveX Client Control must be installed and active.
- The client PC does not need to be a Windows 7 or Windows 2008 Server. Rather, it can be a Windows XP PC, for example.
- If you are using Windows XP for the client PC, be sure to install SP3.




Setting up RD Web Access for IM

Overview of Access Methods

Your users can access IM using two available remote desktop technology methods. Remote desktop technology setup varies slightly, depending upon your chosen access method.

Overview of Windows Installation access

Using a Windows Installation, users can double-click on a desktop or start menu icon to access IM. This access method simulates running IM on the user's PC; however, IM is actually running on the server.

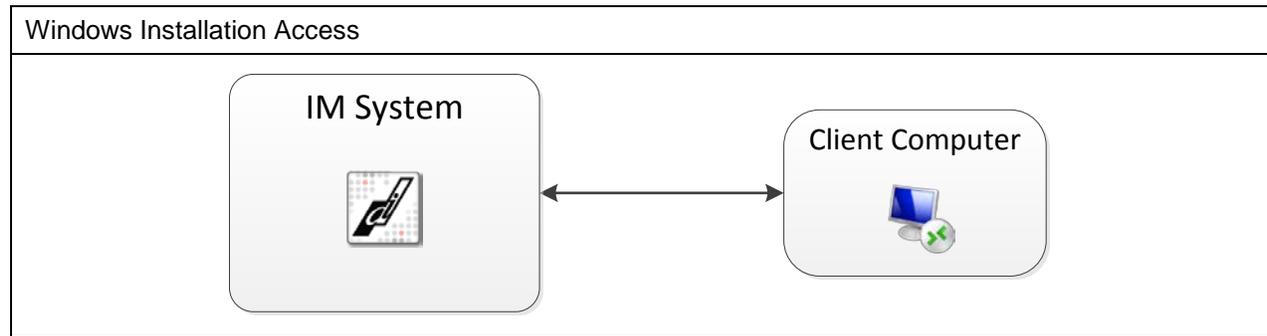


Steps for setting up this access method include:

#	Step / Description
1	Set up the server.
2	On the server, create a Windows Install Microsoft Installer (.msi) file.
3	Install the .msi file on the client PC.

Overview of Web URL access

Using a Web URL, users can access IM through an ActiveX control running inside Internet Explorer's browser to launch Instrument Manager. In this mode, IM still runs on the server.

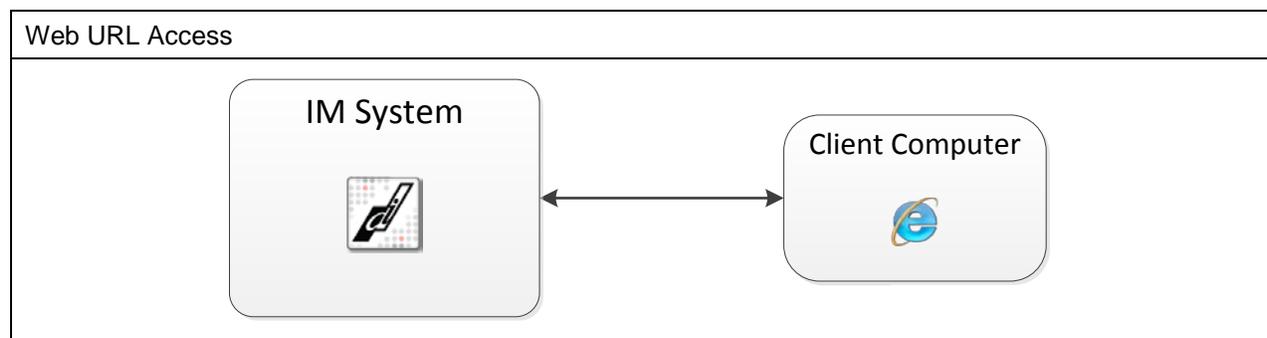

Steps for setting up this access method include:

#	Step / Description
1	Set up the server.
2	On the server, create and modify a Remote Desktop Protocol (.rdp) file.
3	Enable the client PC to connect to the server.



Setting up the Server

To install RD Web Access:

1. On the server, download the **Remote Server Management Toolkit** from Microsoft:
<http://www.microsoft.com/download/en/details.aspx?displaylang=en&id=7887>
2. To install the toolkit, double-click the downloaded file.
3. In the installation wizard dialogue boxes, follow the instructions to install remote desktop tools.

To enable RD Web Access:

To set up RD Web Access on the same server, enable Remote Desktop Services:

#	Step / Description
1	On the server, go to Server Manager > Roles > Add Roles .
2	Click Next .
3	Check Remote Desktop Services and Web Server (IIS) .
4	Click Next .
5	Click Next .
6	Check Remote Desktop Session Host and Remote Desktop Web Access .
NOTE: If you are prompted for required role services, accept the call by clicking Add Required Role Services.	
7	Click Next .
8	Check Do not require Network Level Authentication .
9	Click Next .
10	Check Configure Later .
11	Click Next .
12	Accept that the only user or user groups configured is Administrators .
13	Click Next .
14	On the Configure Client Experience dialogue, accept the defaults and then click Next .
15	On the Web Server (IIS) tab, click Next .
16	On the Roles Services tab, click Next .
17	On the Confirm Installation Selections tab, click Install .
18	Wait for the installation to complete.
19	Click Close .
20	Follow the instructions to shut down the server, then reboot the server.
21	In the final installation completion dialogue box, click Close .
22	Go to Start > Run .
23	In the command prompt, type services.msc , then press Enter .



Set up the Windows Install:

To create a Windows Install .msi file:

#	Step / Description
1	On the server, go to Server Manager > Roles > Remote Desktop Services .
2	Click RemoteApp Manager .
3	In the RemoteApp Manager , in the bottom of the horizontal middle grid, right-click Add RemoteApp Programs .
4	In the RemoteApp Wizard , select Instrument Manager , then click Next .
5	On the next screen, click Finish .
6	In the grid that is displayed, right-click the Instrument Manager entry that you just created, then select Create Windows Installer Package .
7	In the initial Create Windows Installer Package Wizard screen, click Next .
8	In the next wizard dialogue box, do one of the following: <ul style="list-style-type: none"> Note the default location where the wizard saves the installer file (e.g., C:\Program Files\Packaged Programs). Browse to a preferred location where the wizard saves the installer file.
9	Click Next .
10	In the wizard dialogue box, check Desktop and Start Menu Folder .
11	Click Next .
12	Click Finish . A Windows Explorer window is opened to the folder you chose in step 5.

To install the Windows Install .msi file on the client PC:

#	Step / Description
1	From the server folder where you created the Windows Install .msi file, copy the Instrument Manager.msi file from this server.
TIP: For example, to copy and paste the .msi file, use the network or a thumb drive.	
2	Paste the .msi file to the PC that you want to use a client PC.
3	On the client PC, double-click the .msi file to run the Windows installation.
4	Once the installation is complete, double-click the IM icon that the installation automatically put on the desktop. (This is the IM executable file.)
5	In the RemoteApp dialogue box, click Connect .
6	Log in using credentials for the server.

NOTE: If you are prompted that there is a missing certificate (i.e., that the server cannot be identified), ignore the message and click OK.



To verify the Windows Install setup:

#	Step / Description
1	On the client PC, access the IM executable file.
2	Log in to IM as Administrator (IM_ADMIN).
3	Go to System > Status of Instrument Manager Sessions . In the Status of Instrument Manager Sessions screen, note the PC Name and IP Address .
4	On the client PC, go to Start > Run .
5	Type cmd .
6	In the command prompt, type ipconfig , then press Enter .
7	Verify that the PC client name and IP address match those in the session of IM.

Set up the Web URL

To create an .rdp file:

#	Step / Description
1	On the server, go to Server Manager > Roles > Remote Desktop Services .
2	Click RemoteApp Manager .
3	In the RemoteApp Manager , in the bottom of the horizontal middle grid, right-click Add RemoteApp Programs .
4	In the RemoteApp Wizard , select Instrument Manager .
5	Click Next .
6	Click Finish .
7	In the grid that is displayed, right-click the Instrument Manager entry that you just created, then click Create .rdp File .
8	Click Next .
9	In the initial Remote App Wizard screen, click Next .
10	In Specify Package Settings , in Enter the location to save the packages , do one of the following: <ul style="list-style-type: none"> Note the default location where the .rdp file is saved (e.g., C:\Program Files\Packaged Programs) Browse to a new location where you want to save the .rdp file.
11	Click Next
12	In Remote Desktop Session Host Settings , click Change to modify the server name or the Remote Desktop Protocol (RDP) port number .
13	Click OK
14	In RD Gateway Settings , click Change to modify or to configure whether clients will use an RD Gateway server to connect to the target RD Session Host server across a firewall .
15	Click OK
16	In the Certificate Settings , click Change to select or to change the certificate to use .
17	Select the certificate that you want use to digitally sign the .rdp file, and then click OK .



To enable the client PC to connect to the server:

#	Step / Description
1	On the client PC, use Internet Explorer to access: <a href="https://<server-name>/rdweb">https://<server-name>/rdweb
NOTE: In the above URL, replace <server-name> with the hostname of the server where the Instrument Manager application was placed.	
2	If you are prompted for your network login and password, enter that information.
3	Click OK .
4	In the Remote Desktop Services Default Connection screen, do the following: <ul style="list-style-type: none"> Note whether your ActiveX control is installed or not. If you are running Windows XP, follow the instructions to ensure SP3 is installed before continuing.
5	Activate the ActiveX control by the following sequence:
6	In Internet Explorer, go to Tools > Manage Add-Ons > Enable or Disable Add-Ons .
7	In the Disabled list, check Microsoft Terminal Services Client Control (redist) .
8	Select Enable .
9	Click OK .
10	Close Internet Explorer.

To verify the Web URL setup:

#	Step / Description
1	On the client PC, use Internet Explorer to access: <a href="https://<server-name>/rdweb">https://<server-name>/rdweb
NOTE: In the above URL, replace <server-name> with the hostname of the server where the Instrument Manager application was placed.	
2	When prompted, log in. An IM icon is displayed on the website.
3	Click the IM icon. A connection dialogue box is displayed
4	Click Connect .
5	Log in with your network credentials.
6	If the certificate on the remote computer cannot not be authenticated when prompted to continue, ignore this warning and then click Yes .
NOTE: You are now logged into an instance of IM, which is running through the ActiveX control that was already part of Windows	



Accessing IM with RD Web Access

Windows Install

To access IM with RD Web Access via a Windows Install:

#	Step / Description
1	On the client PC, access the IM executable shortcut file.
2	Log in to IM.

Web URL

To access IM with RD Web Access via a Web URL:

#	Step / Description
1	On the client PC, use Internet Explorer to access: <a href="https://<server-name>/rdweb">https://<server-name>/rdweb
NOTE: In the above URL, replace <server-name> with the hostname of the server where the Instrument Manager application was placed.	
2	When prompted, log in. An IM icon is displayed on the website.
3	Click the IM icon. A connection dialogue box is displayed.
4	Click Connect .
5	Log in with your network credentials.